

Redditch Medical InSpec™ DE

Product Description

InSpec™ DE is a sterile preparation of 70% Denatured Ethanol (DE) in 30% Water For Injection (WFI) quality water.

General Guidance

- Use Instructions: InSpec™ DE is designed for spraying, wiping, and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate bucket for mopping. Apply to surface to ensure complete coverage. For 5L screw cap bottles use the entire contents after opening.
- Appearance: Clear liquid free from particles. Specific Gravity (20°C): 0.880 to 0.888. Denatured Ethanol proportion: 70% +/- 2% v/v. This data is typical of normal production and should not be taken as a specification.
- Storage: Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is provided in a separate Safety Data Sheet (SDS).
- Material Compatibility: Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical

Benefits

- Low residue.
- Rapid action.
- Bactericidal and Fungicidal.
- 1 Litre Trigger Spray has 3 months validated 'in-use' shelf life as contents are protected via 'bag-in-bottle' technology.
- All formats manufactured with WFI quality water.
- Multiple bags for cleanroom transfer.
- Ideal surface and aseptic transfer disinfectants.
- Manufactured in accordance with GMP.



Technical User Information

- InSpec™ DE is a sterile ready to use disinfectant. This is a 70% v/v denatured ethanol (TSDA 7) solution in WFI quality water.
- InSpec™ DE is manufactured in accordance with GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.
- The solution is filtered aseptically through a 0.2-micron filter and gamma irradiated to a validated process (at 25kGy) to give a sterility assurance level SAL of 10⁻⁶.
- InSpec™ DE comes with key certification such as irradiation, sterility, endotoxin, and analysis.

Purchase Information

DEWFIAUS-LS

DEWFIAUS-500MLS

Packaging

6 x 1 L Trigger spray bottles/case

8 x 500 mL Trigger spray bottles/case

