

Redditch Medical InSpec™ QT

Product Description

High performing broad-spectrum disinfectant for hard surfaces. QT+ is a low residue twin chain quaternary ammonium compound with cleaning and disinfecting properties.

General Guidance

- Use Instructions: InSpec™ QT is designed for spraying, wiping, and mopping applications. Hold spray approximately 15cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. 100ml concentrate added to 2.5L of water is required for mopping. Apply to surface to ensure complete coverage.
- Technical Data: Aldehyde-free disinfectant cleaner. 4% RTU Solution pH 11.8-12.2 Concentrate pH 12.9. This data is typical of normal production and should not be taken as a specification.
- Storage: Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is provided in a separate Safety Data Sheet (SDS).
- Material Compatibility: Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical Dossier.

Benefits

- An exceptional non-oxidising biocide.
- Bactericidal and fungicidal.
- Clean and disinfects in the presence of dirt, blood, and proteins.
- 3 months validated 'in-use' shelf life.
- Contents protected using 'bag-in-bottle' presentation with premium range.
- 1L and 5L formats manufactured with WFI quality water.
- Multiple bags for cleanroom transfer.
- Ideal rotational partner for InSpec AN, HA and OX.
- Manufactured in accordance with GMP.



Technical User Information

- InSpec™ QT is a sterile disinfectant in pre-diluted and concentrated formats. InSpec™ QT is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.
- The solution is filtered through a 0.2-micron filter and gamma irradiated to a validated process (at 25-45kGy) to give a sterility assurance level SAL of 10⁻⁶.
- Trigger sprays supplied with 900ml solution in a 1L container.
- InSpec™ QT is supplied with Certificates of Conformity, Analysis, Irradiation Sterility and Endotoxin.
- All formats of QT have a stabilised 24-month shelf life.
- For screw cap bottles use the entire contents after opening.

Purchase Information

QTPLWFIAUS-900MLS

QTPLCNCAUS-102

Packaging

6 x 900 mL Trigger spray bottles/case

5 x 100 mL bottles per double bagged pack,
10 packs per case, total 50 bottles per case



MMRC Pty Ltd
Cleanroom & Contamination Control

ABN: 41 650 603 295

P. +61 3 9885 1752
E. mail@mmrc.com.au

www.mmrc.com.au

11 Thornton Cres, Mitcham
VIC 3132 Australia