



Technical Datasheet

AN



Sterile Diamine Disinfectant Cleaner

Low residue QUAT free disinfectant / cleaner. Broad spectrum disinfectant.

Use Instructions

InSpec™ AN is designed for spraying, wiping and mopping applications. Hold spray approximately 1.5cm to 20cm from area to be treated. 5L versions, pour into an appropriate container for mopping. 100ml concentrate added to 2.4L of water is required for mopping. Apply to surface to ensure complete coverage.

Technical Data

Alcohol, chlorine and QUAT-free disinfectant cleaner. 4% RTU Solution pH 8.8-9.2, Concentrate pH 10.0.

Storage

Store upright in original closed containers, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is available in the Safety Data Sheet (SDS).

Safe Handling

Always wear gloves and goggles or face protection. Always read the label and SDS before use.

Material Compatibility

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.

Benefits

- An exceptional non-oxidising biocide
- Bactericidal and Fungicidal
- Clean and disinfects
- 3 months validated 'in-use' shelf life
- Contents protected using 'bag-in-bottle' presentation with premium range.
- 1L and 5L formats manufactured with WFI quality water
- Multiple bags for cleanroom transfer
- Ideal rotational partner for **InSpec™** QT, HA and OX
- Manufactured in accordance with GMP

Technical User Information

InSpec™ AN is a sterile disinfectant in pre-diluted and concentrated formats.

InSpec™ AN is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter and gamma irradiated to a validated process (at 25-45kGy) to give a sterility assurance level SAL of 10⁻⁶.

Trigger sprays supplied with 900ml solution in a 1L container.

InSpec™ AN is supplied with Certificates of Conformity, Analysis, Irradiation and Sterility.

All formats of **InSpec™ AN** have a stabilised 24 month shelf life. For screw cap bottles use the entire contents after opening.

Microbiological Minimum Efficacy Testing Contact

EN 1276 Bacteria	5 minutes	EN 13697 Bacteria	5 minutes
EN 1650 Fungi	15 minutes	EN 13697 Fungi	15 minutes

InSpec™ AN Product

Product	Pack Size	Case Size	Product Code
InSpec™ AN 900mls Premium - Sterile	900mls trigger spray	6 x 900mls	ANWFI20-1LS
InSpec™ AN 5L Screw Cap - Sterile	5L Screw cap	2 x 5L	ANWFI20-5LS
InSpec™ AN Concentrate - Sterile	100ml Concentrate	50 x 100mls	ANCNC20-100
InSpec™ AN Bustable Pouch -Sterile	15 Wipes/pouch	20 x 15	ANWFI20BP15S
InSpec™ AN 900mls Premium - Non-Sterile	900mls trigger spray	6 x 900mls	NSANWFI-1L
InSpec™ AN 5L Screw Cap - Non-Sterile	5L Screw cap	2 x 5L	NSANWFI-5L
InSpec™ AN Concentrate - Non-Sterile	100ml Concentrate	50 x 100mls	NSANCONC-100ML